


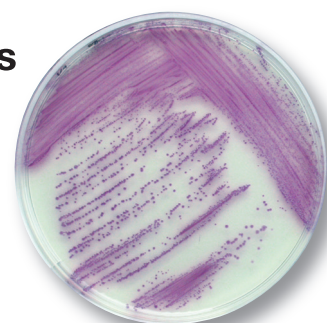

RAPID'*Salmonella* Short Protocol

- **Detection of *Salmonella* spp. in less than 42 hours**
- **Easy protocol**
1 single broth, 1 single solid agar plate
- **Easy confirmation**
Salmonella Latex
- **Meets the regulations**
Detects *Salmonella* typhi/paratyphi, lactose-positive, motile and non-motile *Salmonella*
- **Validated medium that can be used as a second medium in the ISO 6579 Standard**




The Difference is Clear. RAPID'Chromogenic Range for Food and Environmental Testing

BIO-RAD

RAPID'Salmonella Short Protocol

Principle

The principle behind RAPID'Salmonella relies on the demonstration of two enzymatic activities. *Salmonella* spp. take the form of typical, readily identified magenta colonies (detection of C8 esterase). Counter selection is used to reveal other bacteria with a different color.

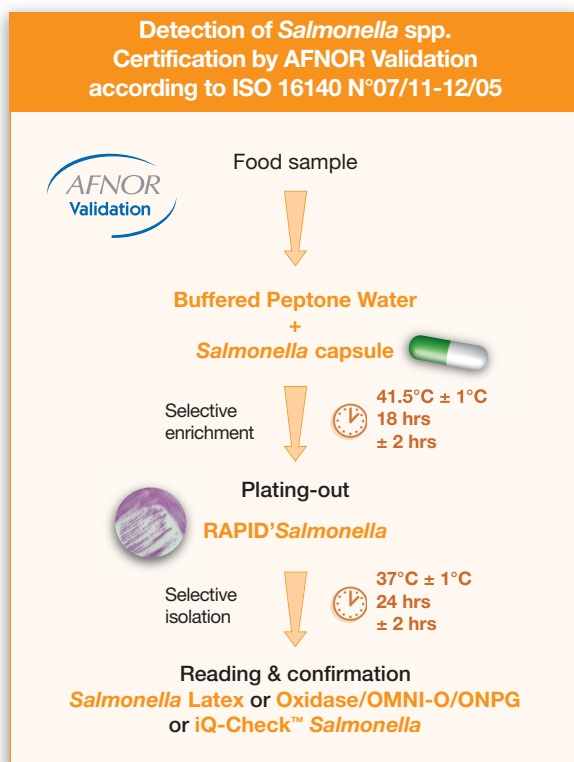

Ordering information

Catalog #	Description
RAPID'Salmonella/Agar	
Pre-poured	
356-3961	20 x 90 mm
Dehydrated	
356-4705	500 g
Buffered Peptone Water	
Dehydrated	
356-4684	500 g
Ready to use	
355-4179	225 ml x 6 bottles
355-5789	2.3 l x 5 bags
355-5790	5 l x 2 bags
Enrichment Supplement	
356-4710	Salmonella capsule100
Confirmation	
355-3832	Oxidase discs30
356-0781	Omni-O (A-60)3 ml
355-3822	ONPG50 discs
355-6710	Latex Salmonella kit75

Key Benefits

Fast results

- Less than 42 hrs for results to detect *Salmonella* spp.
- Fast confirmation (less than 1 minute)

One broth + one plate

- Saves money
- Saves handling time

BPW enrichment can be used for sample dilutions

- No need for another stock solution (such as tryptone salt)

Easy to use

- Solid agar plates
- Easy confirmation

Very low false positive rate (<0.3% all matrices*)

- Saves on useless confirmations

Meets regulatory expectations

- RAPID'Salmonella detects motile and non-motile *Salmonella*, as well as lactose-positive *Salmonella*, *Salmonella typhi* and *Salmonella paratyphi*.

* during ISO 16140 accuracy tests

BIO-RAD

**Bio-Rad
Laboratories, Inc.**

Life Science
Group

Web site www.bio-rad.com **USA** 800 4BIORAD **Australia** 61 02 9914 2800 **Austria** 43 (0) 1 877 89 01 **Belgium** 09 385 55 11 **Brazil** 55 21 3237 9400
Canada 1 800 268 0213 **China** 86 21 6305 2255 **Czech Republic** 420 241 430 532 **Denmark** 44 52 10 00 **Finland** 09 804 22 00 **France** 33 1 47 95 62 59
Germany 49 (0) 89 318 84 0 **Greece** 30 210 777 4396 **Hong Kong** 852 2789 3300 **Hungary** 36 1 455 8800 **India** 91 124 2398 112/113/114
Israel 03 951 4127 **Italy** 39 02 216 091 **Japan** 03 5811 6270 **Korea** 82 2 3473 4460 **Latin America** 305 894 5960 **Mexico** 52 555 200 0520
The Netherlands 31 318 540 666 **New Zealand** 64 9 415 2280 **Norway** 23 38 41 30 **Poland** 48 22 331 99 99 **Portugal** 351 21 472 7700
Romania 4021 210 1703 **Russia** 7 095 721 1404 **Singapore** 65 6415 3188 **South Africa** 27 11 442 8508 **Spain** 34 91 590 5200 **Sweden** 46 8 555 12700
Switzerland 41 (0) 61 717 95 55 **Taiwan** 886 2 2578 7189 **Thailand** 662 651 8311 **United Kingdom** 44 20 8328 2000 **Vietnam** 848 823 6757